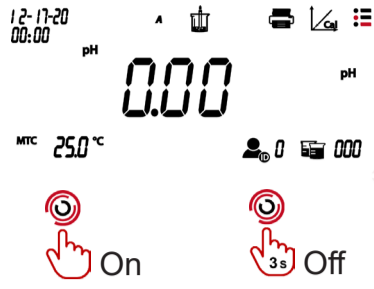
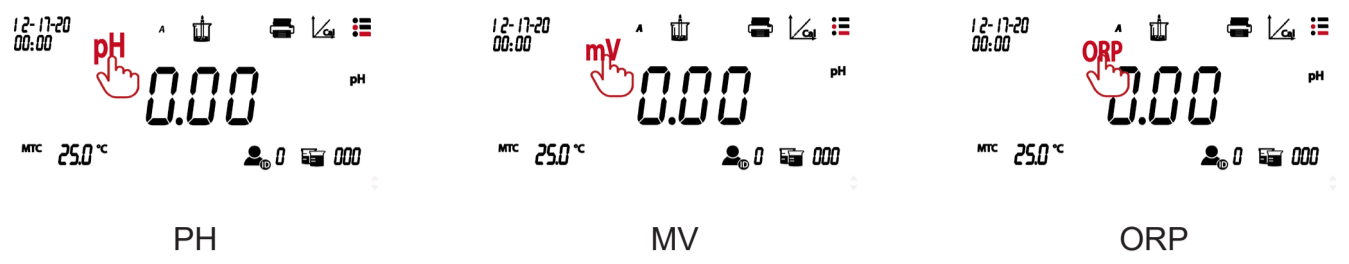
Switch On/Off

Switch pH/mV/ORP Mode

Temperature Unit Setup

Temperature Setup

pH Buffer Solution Setup

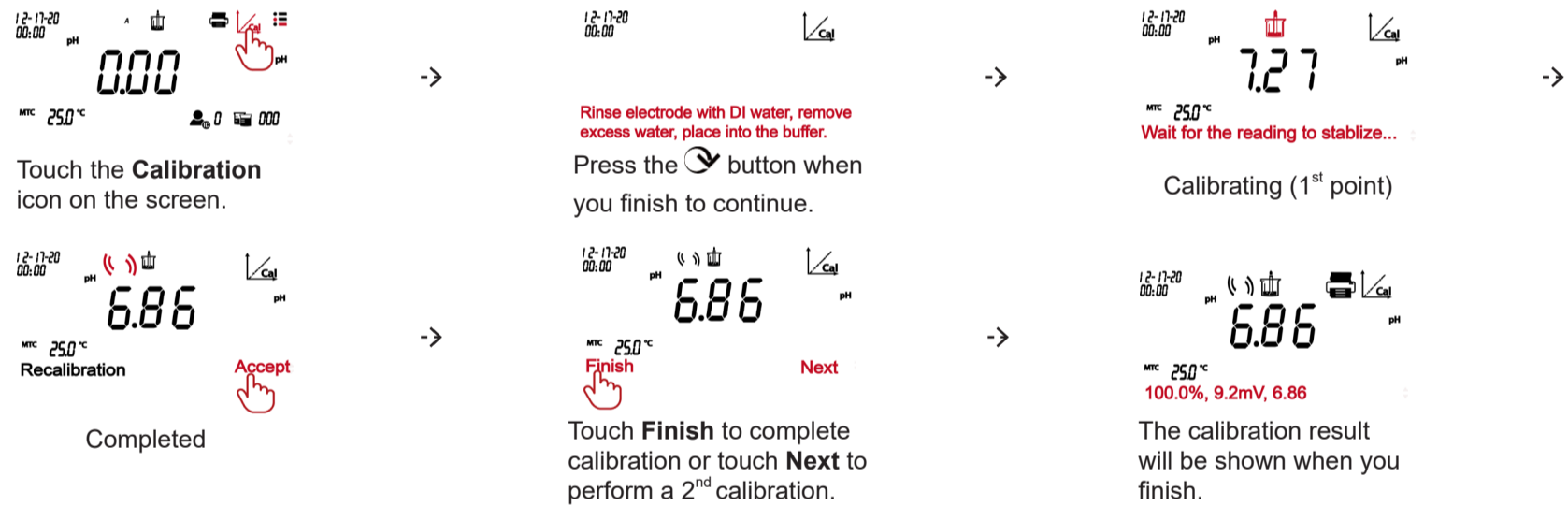
Note: Please check the pH buffer group table below for details of the buffer values.

pH Resolution Setup


Note: pH Resolution value varies for a-AB41PH and a-AB33PH, so please set it based on the actual value on the screen.

Calibration


Measurement

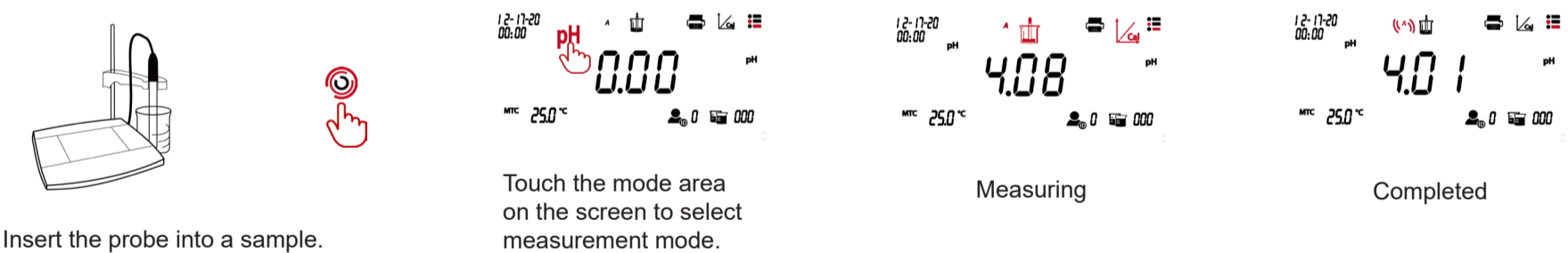
pH Buffer Group

pH Buffer Group	Temperature	pH Value				
USA standard	25°C	1.68	4.01	7	10.01	/
J.J.G standard	25°C	1.68	4.01	6.86	9.18	12.46
Europe standard	25°C	2	4.01	7	9.21	11
MERCK (Ref.20°C)	20°C	2	4	7	9	12
DIN (19266)	25°C	1.68	4.008	6.865	9.184	12.454

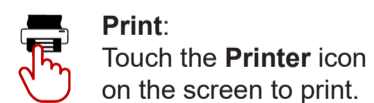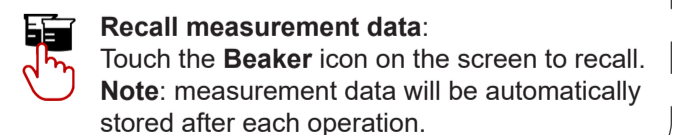
Print


Store/Recall Data

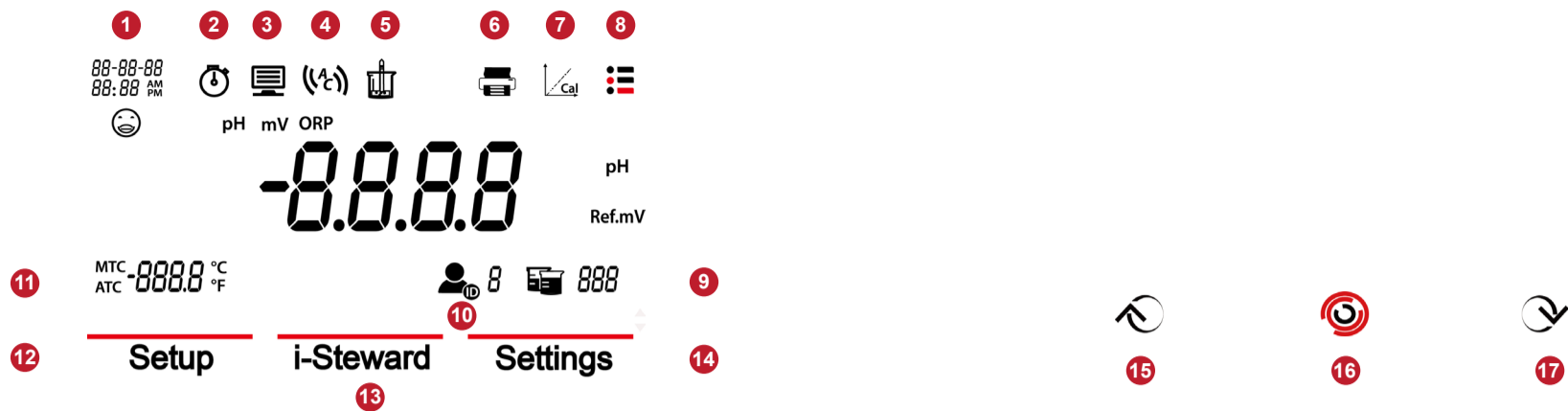Display Introduction

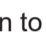1. **Date and time**
2. **Clock icon:** appears when the timing function is used.
Note: this function is only available for a-AB41PH model.
3. **Computer icon:** appears when connects to a computer.
4. **Endpoint mode**
5. **Measurement icon:** blinks when measurement is in progress.
6. **Printer icon:** touch to print.
7. **Calibration icon:** touch to calibrate; blinks when calibration is in progress.
8. **Menu icon:** touch to enter menu settings.
9. **Beaker icon:** touch to recall measurement data.
10. **User ID icon:** touch and then press  or  button to select user by user ID.
11. **Temperature:** touch to switch between MTC and ATC temperature.

Touch the **Menu icon**  to enter the following setup menu:

12. **Setup:** touch to enter parameter setup menu.
13. **i-Steward:** touch to enter sensor condition menu.
14. **Settings:** touch to enter meter's setup menu.

Note: when you enter the three menus introduced above, touch the related sub-menu items to finish settings. Please refer to the first page of the flyer for detailed information about some of the sub-menu items.

Buttons:

15. Up button:

- a. When setting parameters, increases the setting value or displays the previous option.
- b. When in setup menu, press to scroll up to the previous page of menu selections.
- c. When in calibration and PH ECS process, press to repeat the previous step.

16. Home button:


- a. Press to turn on the meter.
- b. Press to start a new measurement.
- c. Press to exist from the current mode or menu and return to the measurement mode.

17. Down button:


- a. when setting parameters, decreases the setting value or displays the next option.
- b. when in setup menu, press to scroll down to the next page of menu selections.
- c. When in calibration and PH ECS process, press to move to the next step.

Install the stand-alone electrode holder

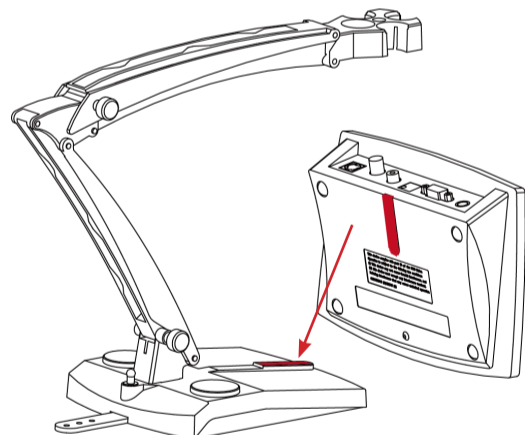
1. Attach the magnetic part on the bottom of the meter to the magnetic part on the stand-alone electrode holder to finish installation. (Both marked in red demonstrated below.)

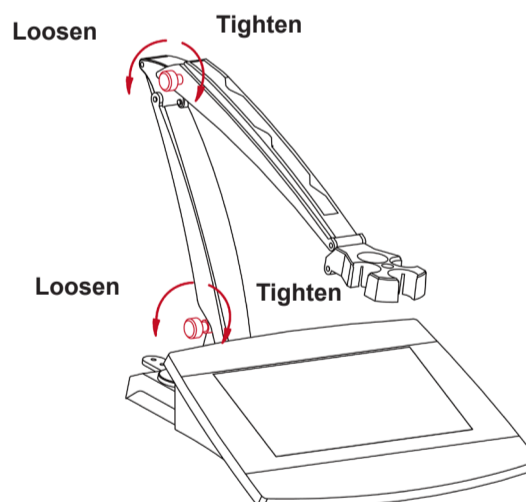2. The two tension knobs can be adjusted by turning clockwise to tighten or turning counterclockwise to loosen. When they are loosened, you can move the holder up and down, forward and backward to the desired angle. Then you can tighten the knobs to fix the holder.

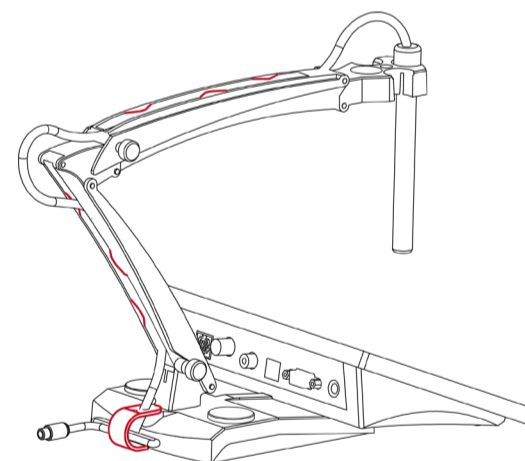
Note: the stand-alone holder boasts a harness storage design. You can push the electrode's cable in the slots marked in red below, roll up and fix the harness storage strip located at the bottom of the holder, and pass the cable through the strip.

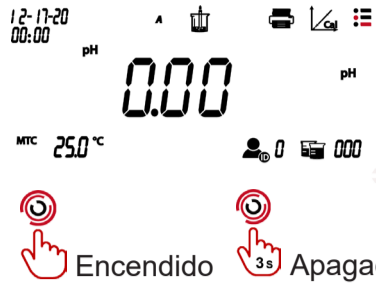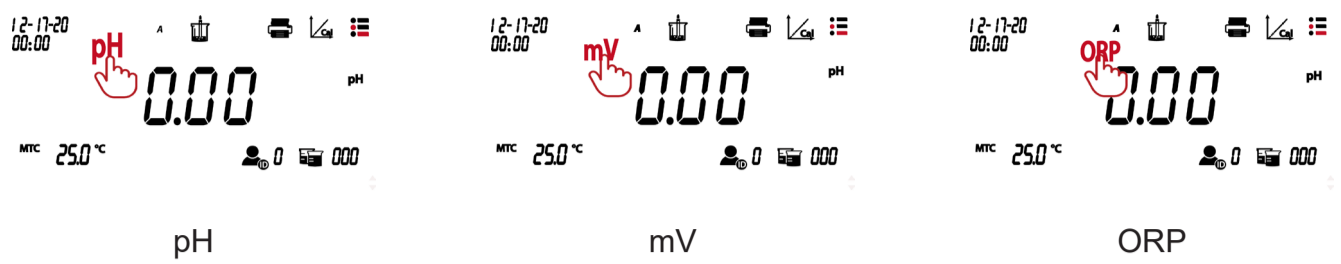
Scan the QR code to get **a-AB33PH** and **a-AB41PHs'** instruction manual.

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