



GUARDIAN™ SERIES

Hot Top Indicator Guide

Think Safe. Think SMART.

The Guardian™ series Hotplates and Stirrers feature a unique safety feature to protect users in the lab. Knowing when the top plate is hot to the touch can prevent accidental burns. To keep users safe, all Guardian models have a hot top indicator light.

This 13 mm hot top indicator light is prominently featured on all models from the Guardian 2000 to the Guardian 7000. Once the Guardian's heater reaches 40°C, the indicator light turns on and will stay illuminated even when the Guardian power switch is turned off. Only when the Guardian's heater temperature drops below 40°C will the indicator light shut off.



Ingeniously Practical

GUARDIAN™ SERIES *Hot Top Indicator Guide*

From the basic Guardian 2000 to the intelligent Guardian 7000, every Guardian has a prominent hot top warning light.

The Guardian series features 4 levels of hotplates and stirrers, the Guardian 2000, 3000, 5000, and 7000.



GUARDIAN 2000



GUARDIAN 3000



GUARDIAN 5000



GUARDIAN 7000

The hot top indicator always illuminates once the heater reaches 40°C

Guardian Series Heated Under 40°C



GUARDIAN 2000



GUARDIAN 3000



GUARDIAN 5000



GUARDIAN 7000

Guardian Series Heated Over 40°C



GUARDIAN 2000



GUARDIAN 3000



GUARDIAN 5000



GUARDIAN 7000

GUARDIAN™ SERIES *Hot Top Indicator Guide*

The hot top indicator will stay lit when over 40°C even when the power is shut off.

Once the power buttons are switched off, the screens for the Guardian 3000, 5000, and 7000 will change to display prominently the word "HOT". Additionally, on the Guardian 5000 and 7000, the screen displays the current heater temperature as the unit is cooling off. The analog Guardian 2000 is without display, yet the hot top indicator light also remains lit after the power is switched off.



GUARDIAN 2000



GUARDIAN 3000



GUARDIAN 5000



GUARDIAN 7000

The warning light stays lit until it cools below 40°C

Once the Guardian reaches a safe temperature, all hot top indicator features will shut off. The Guardian heating surface has now cooled off so is safe if touched by the user. (The Guardian must be plugged in for this feature to work)



GUARDIAN 2000



GUARDIAN 3000



GUARDIAN 5000



GUARDIAN 7000



Click to Watch the Guardian Hot Top Indicator Guide Below



OHAUS Corporation

Headquartered in Parsippany, NJ, OHAUS Corporation manufactures an extensive line of balances and scales, lab equipment and lab instruments that meet the weighing, sample processing and measurement needs of multiple industries. We are a global leader in the laboratory, industrial and education markets, as well as a host of specialty markets, including the food preparation, pharmacy and jewelry industries. An ISO 9001:2015 manufacturer, OHAUS lab balances, industrial scales, lab equipment and lab instruments are precise, reliable and affordable, and backed by industry-leading customer support.

OHAUS CORPORATION
7 Campus Drive
Suite 310
Parsippany, NJ 07054 USA

Tel: 800.672.7722
973.377.9000
Fax: 973.944.7177
www.ohaus.com

The management system governing the manufacture of this product is ISO 9001:2015 certified.



Ingeniously Practical